

Whisper Zoom

Whisper Zoom is a new class of methods designed for highly sensitive analysis of low sample amount.

STEP-BY-STEP GUIDE

- 1 Select the "Instrument configuration" method and run the "Set to Zoom" script: This sets the instrument to the Zoom Loop Mode configuration.
- 2 Please follow the on screen prompt from the script. The transfer line must be connected in port 3 of the Loop valve.
- 3 Please follow the on screen prompt from the script. The HP flow sensor line must be connected in port 6 of the Loop valve.
- 4 Re-tighten the connections.
- 5 Generate and run your sample list using your preferred Whisper Zoom methods.

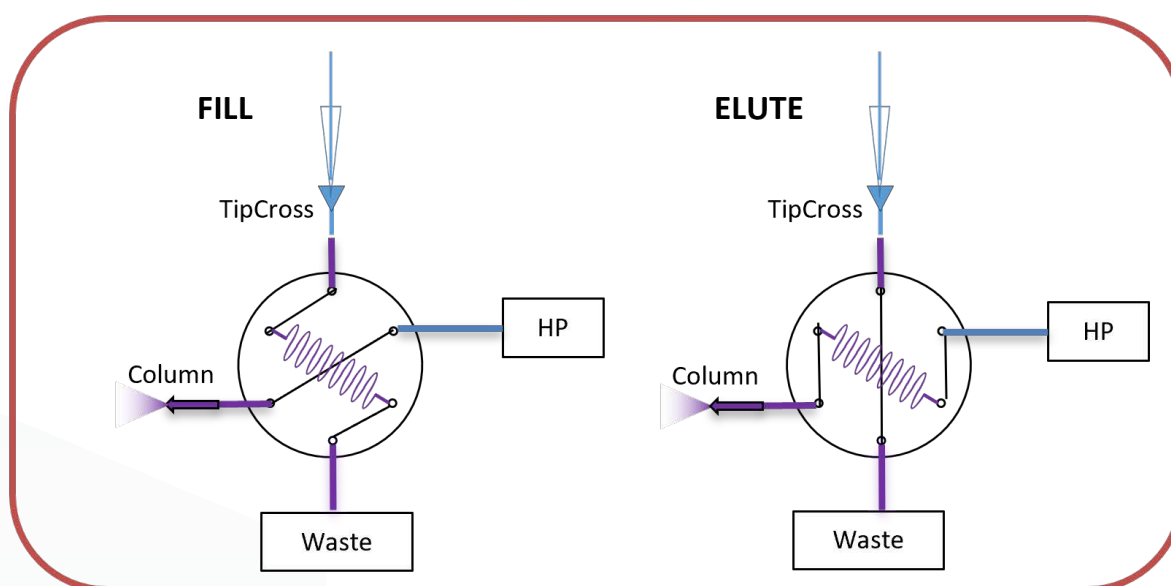


Figure 1: Whisper Zoom configuration